

A COMPACT, HIGH-RESOLUTION TIME-OF-FLIGHT MASS SPECTROMETER BASED ON AN ELECTROSTATIC ANALYSER.

Steven Daly¹, Jan Commandeur¹, Luis Lopez², Kristina Lorenzen², John Hoyes³, Charlotte Uetrecht^{4,5}




1. MS Vision, Televisieweg 40, 1322AM, Netherlands.
2. Scientific Instrument SPB/SFX, European XFEL GmbH, Holzkoppel 4, 22869, Germany.
3. HGSG Ltd., Rowarth, UK.
4. Centre for Structural Systems Biology (CSSB), Deutsches Elektronen-Synchrotron DESY & Leibniz Institute of Virology (LIV), Notkestraße 85, 22607, Germany.
5. University of Siegen, Am Eichenhang 50, 57076, Siegen, Germany




Introduction

Design constraints.

- A diagnostic mass spectrometer was required for the MS-SPIDOC project. Must be able to characterise ion beam containing biomolecular samples including virus capsids.
- Specifications:
 - Fit into small footprint (right).
 - Resolution > 5,000
- Reflectron TOF will not fit → need a different design → ESA.

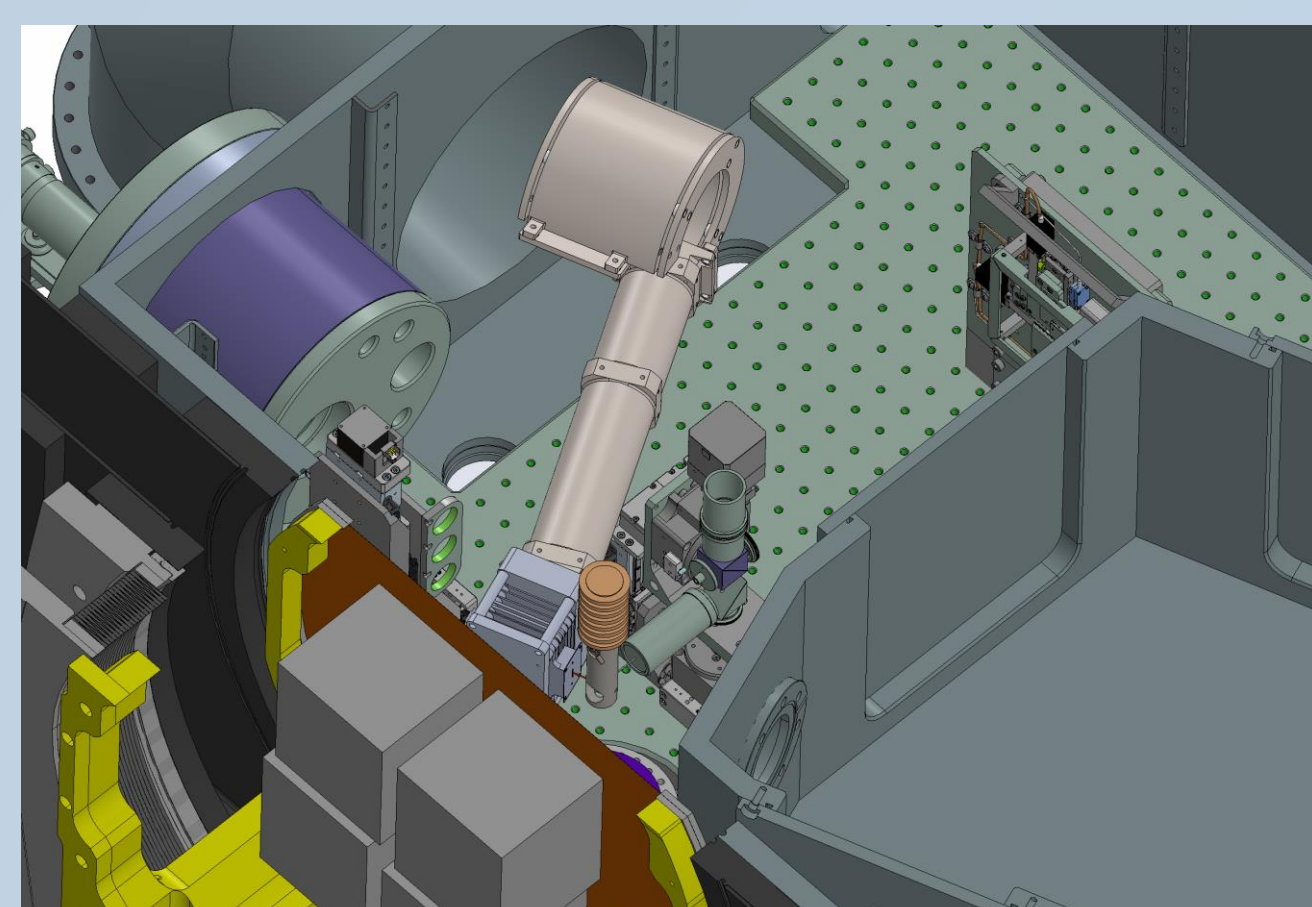

Cutaway view of MS-SPIDOC chamber showing footprint where TOF must fit.

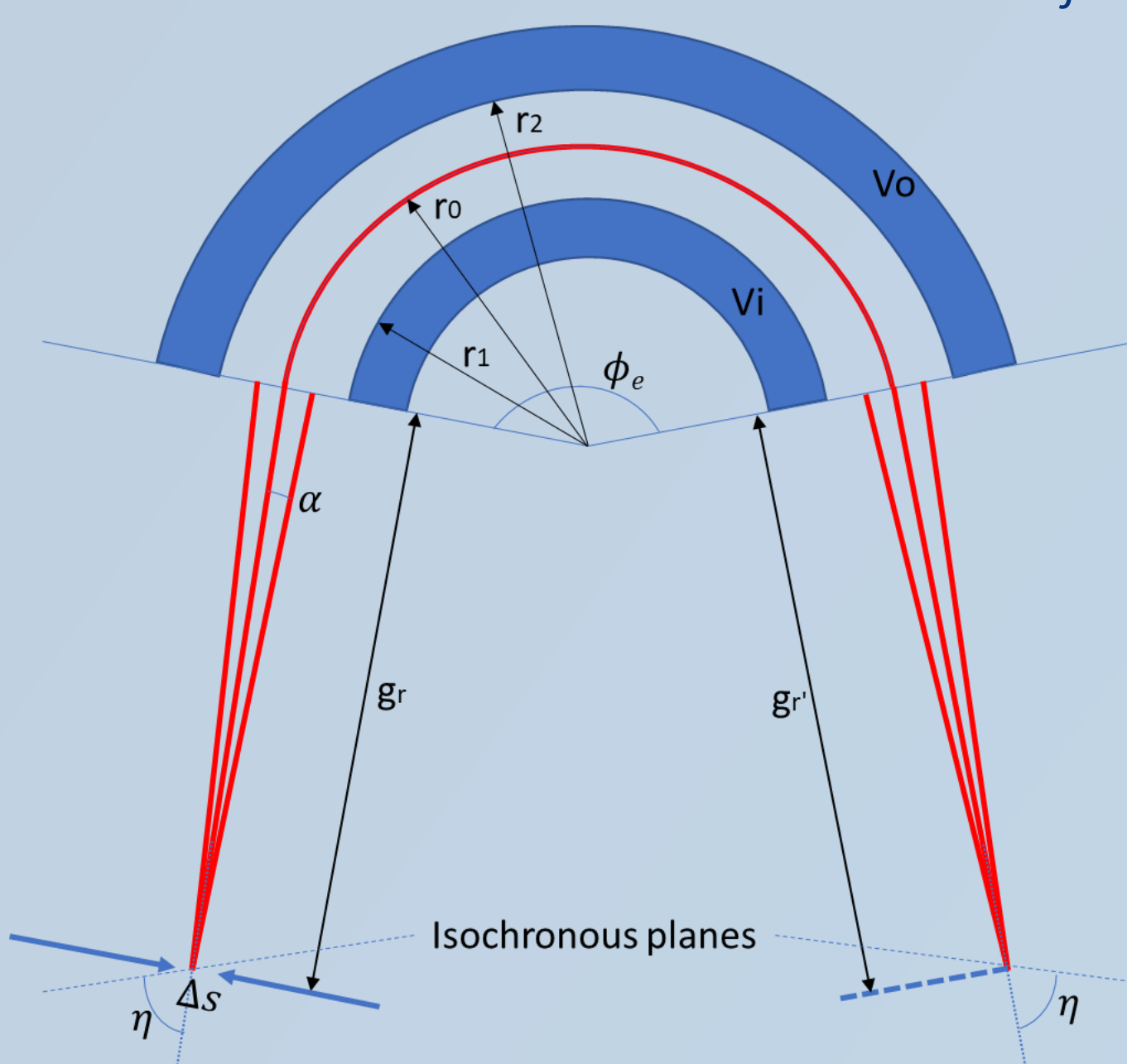


Generalized cylindrical geometry Sector TOF (symmetrical FFR regions).


1. Poschenrider, W.P., Int. J. Mass Spectrom. Ion Processes, 9, 1972, 357-373, [https://doi.org/10.1016/0020-7381\(72\)80020-2](https://doi.org/10.1016/0020-7381(72)80020-2)

- Time of flight mass spectrometers with electrostatic sectors described theoretically in 1972¹.
- When combined with orthogonal acceleration, can arrange the geometry such that second order energy aberration from pusher and ESA cancel.
- Since the ESA acts as an energy filter, it will inherently reject metastable ions, leading to cleaner signal.
- Flexible design → as long as the $g_r + g_r'$ is kept fixed, free to choose the values of g_r and g_r' .

Results: Mass Spectrometry

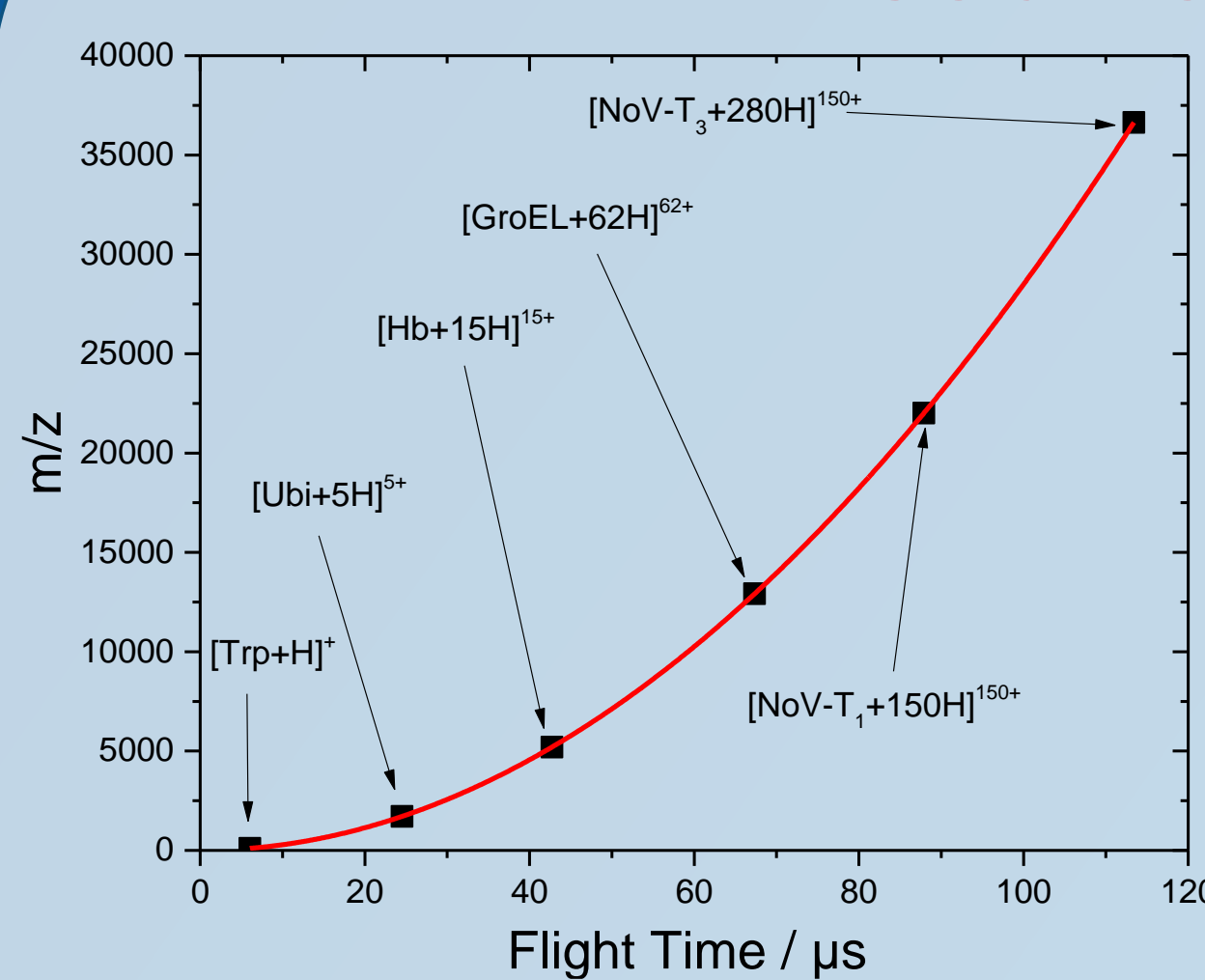


m/z vs flight time for different ions. m/z proportional to TOF squared, as expected.

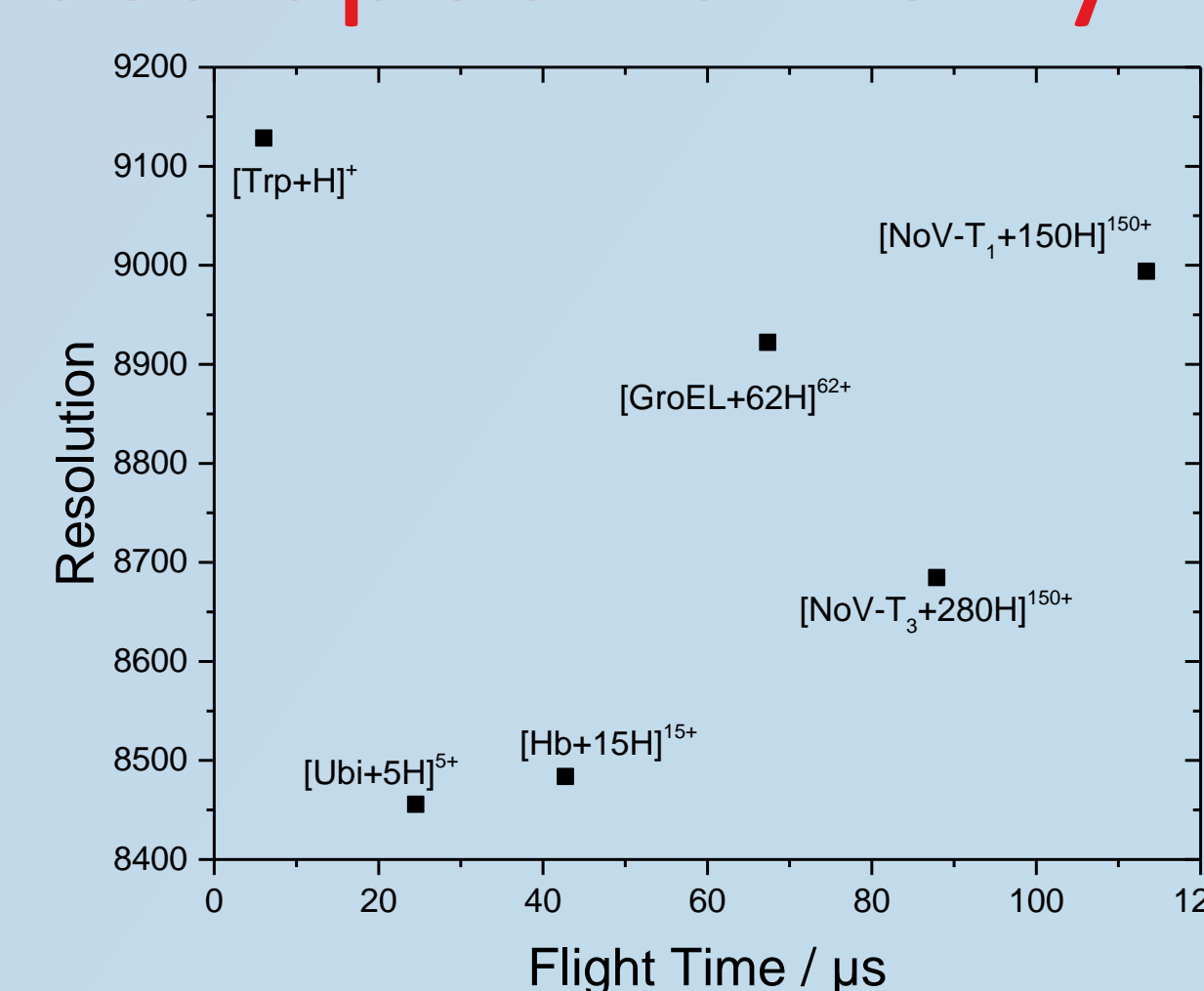


Resolution for ions of different m/z

- Resolution > 8,000 for 23 meV/q energy spread (random).
- For large systems (e.g. viruses), will not be limiting factor in observed resolution.

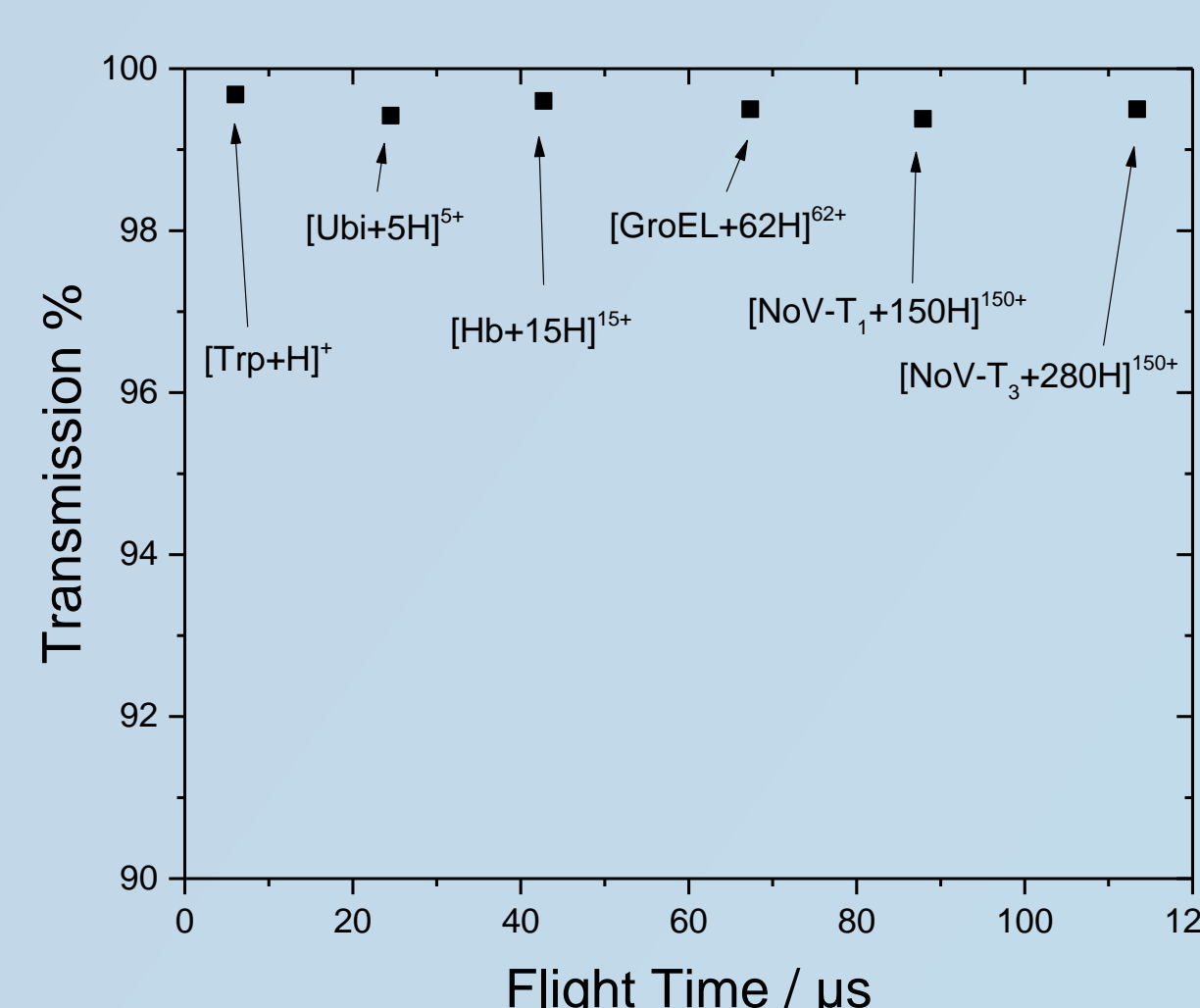


Percentage of ions hitting the detector for different m/z. Transmission > 99% for 23 meV/q energy spread (random). Energy spread in y is main cause for loss of ions on entry to ESA.

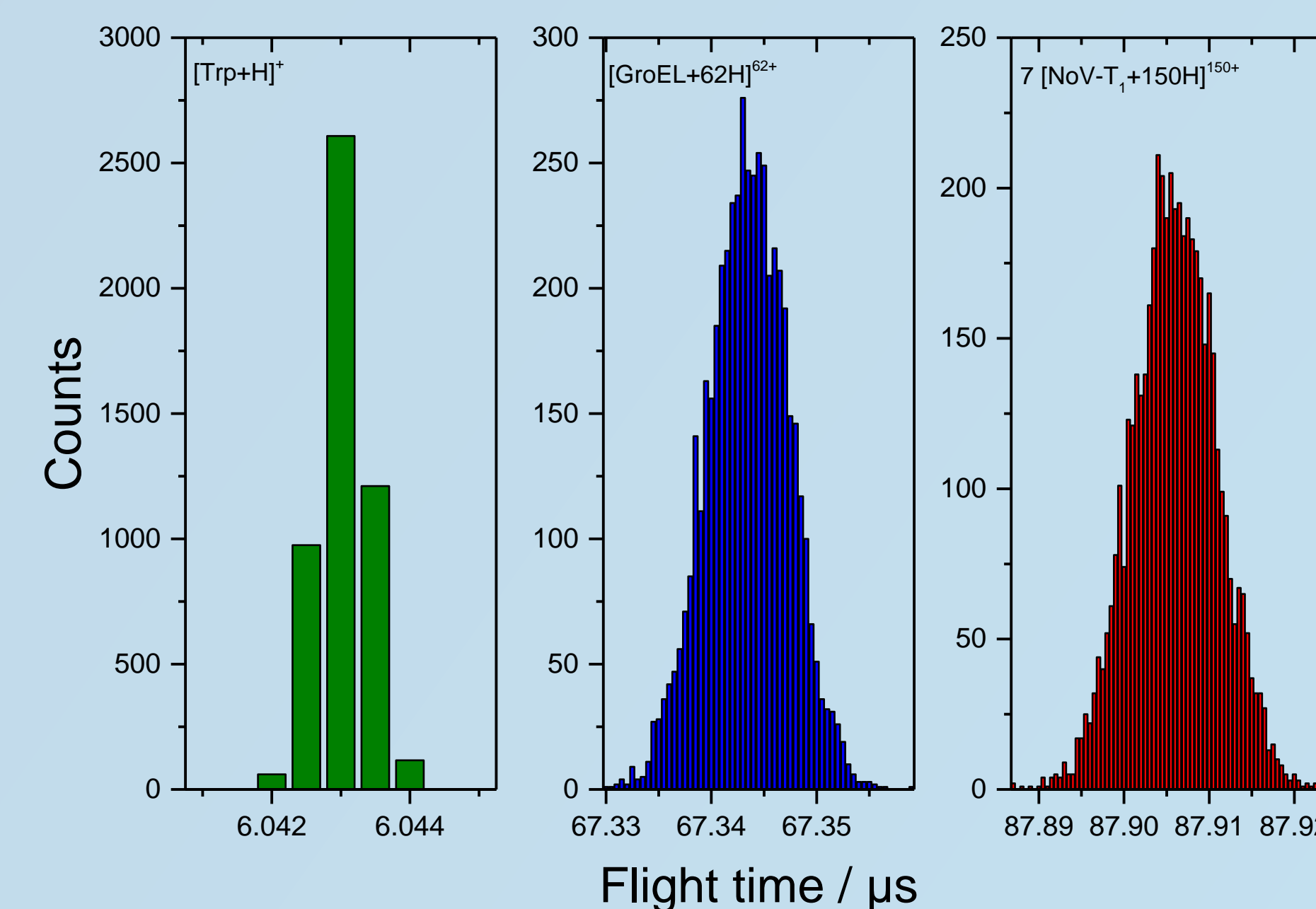

Arrival time distributions at time resolution of ADC for monoisotopic ions with 23 meV/q energy spread (random, 5000 ions).

MS-SPIDOC ESA-TOF Design

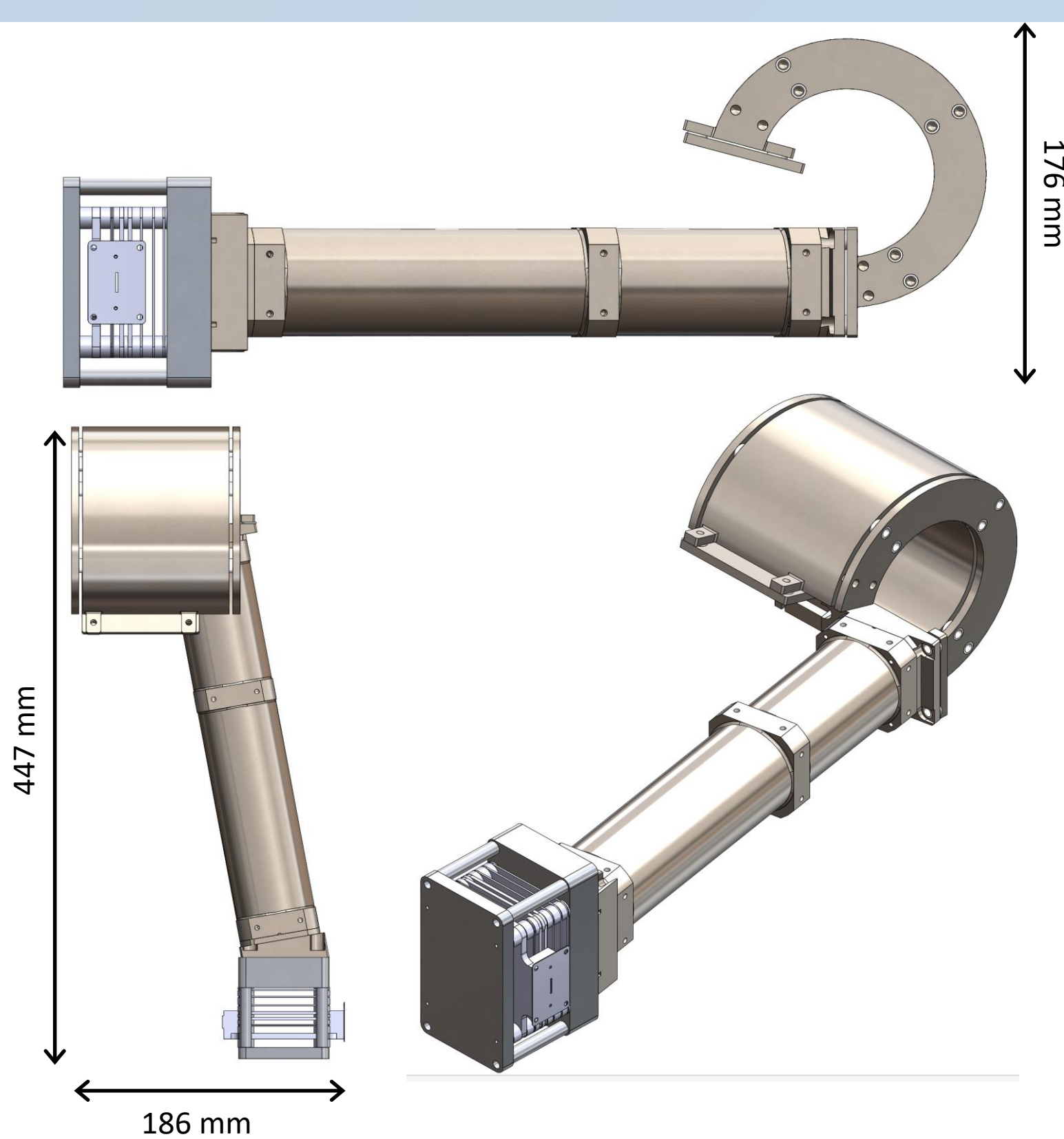

Design of the ESA-TOF for MS-SPIDOC.

Pusher

- Modified QTOF Ultima pusher stack.
- Ions have 1mm spread and 200 eV/q through pusher
- Ions have 6386 eV/q after pusher and travel at 10° to push direction.

Field Free region

- Total length = 244.4 mm, all before ESA in this configuration.


Electrostatic Sector

- $r_1 = 52.25$ mm, $r_2 = 57.75$ mm, $\phi = 254.6^\circ$ (ensure isochronous planes normal to ion trajectories).
- Inner sector has negative voltage relative to FFR, outer sector positive.
- Ion beam traverses 90.4 mm to detector. Ion beam is 25 mm so 27.6 % duty cycle.


Detector

- Ultrafast electron multiplier detector.
- ADC with 500ps resolution


Results: Ultimate Resolution

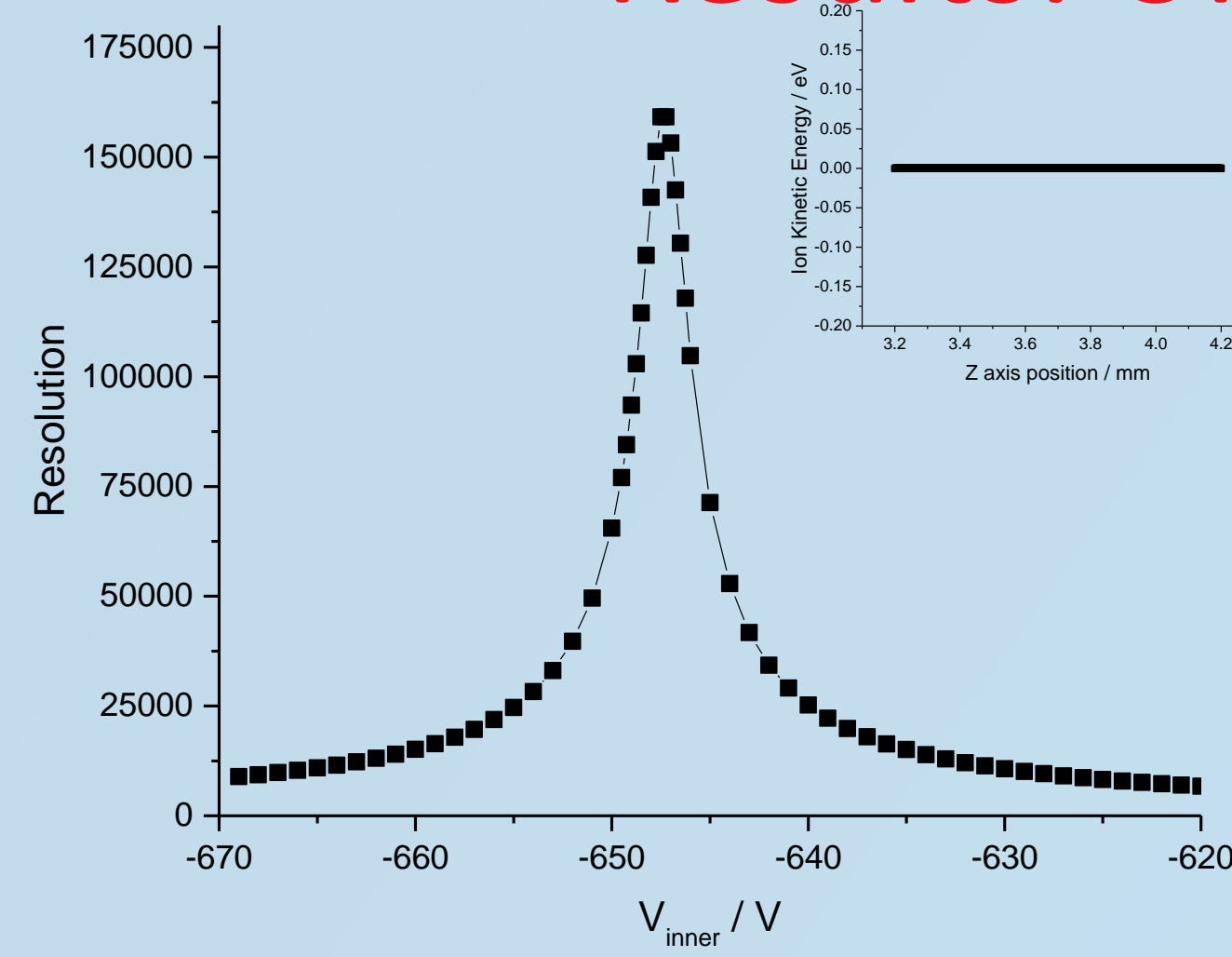

Resolution vs V_{inner} with $V_{outer} = 623$ V for 0 meV/q ion energy spread. Inset shows initial ion distribution in E_z/z plane

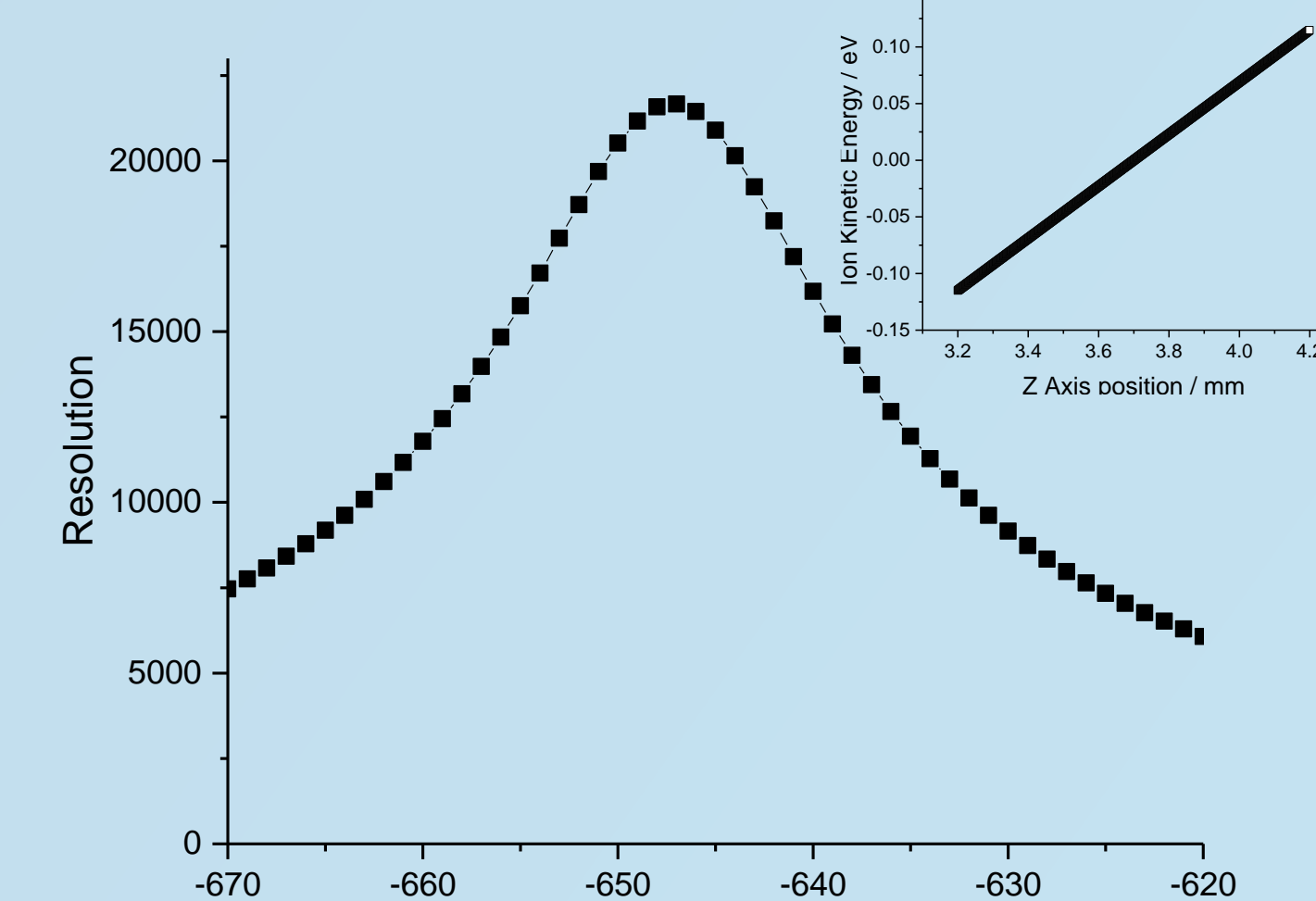

Resolution vs V_{inner} with $V_{outer} = 623$ V for 23 meV/q perfect correlation ion energy spread (inset).

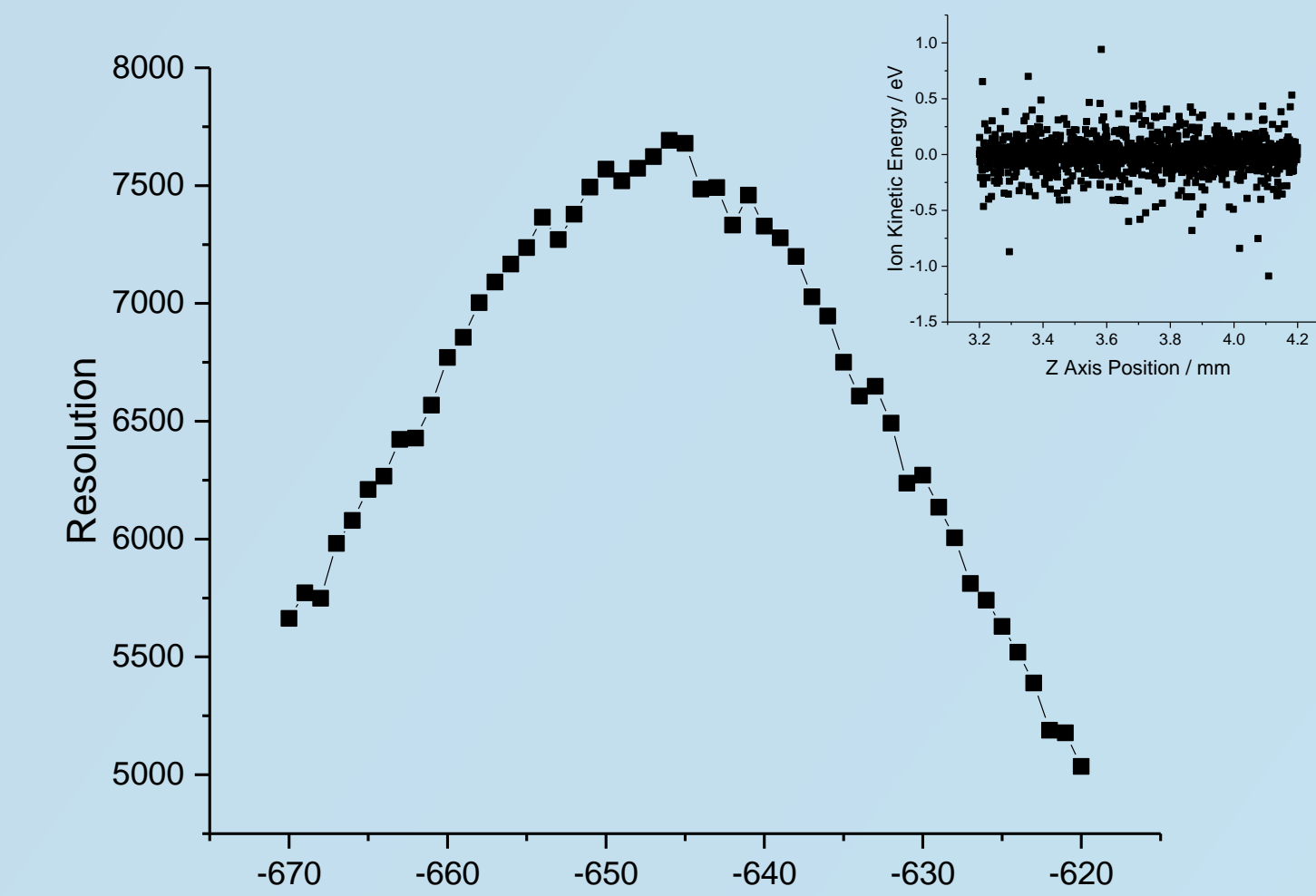

Resolution vs V_{inner} with $V_{outer} = 623$ V for 23 meV/q uncorrelated ion energy spread.

- Very high resolution without any energy spread → detector at second space focus.


- Position and velocity correlated → ions closer to pusher are moving towards & vice versa. Resolution = 22,500 for perfect correlation.

- Fully uncorrelated position and velocity → worst case scenario. Resolution = 7,750.
- Expected performance somewhere between best and worst case → 7,750 – 22,500.

Computational Methods

- Models generated with Simion:
 - 0.05 gu/mm for all.
 - Pusher uses symmetry plane
 - ESA extended 2D model covering entire work bench → no fringe plates needed
 - FFR is empty workbench, only drawn near ESA.
- Simulations performed in SIMAX.


- For all simulations, ions have:
 - $x = 0 \pm 10$ mm, $E_x = 200 \pm 0.023$ eV/q
 - $y = 0 \pm 1$ mm, $E_y = 0 \pm 0.023$ eV/q
- 3 ion distributions in z were used (see resolution).
 - $z = 0 \pm 0.5$ mm, $E_z = 0$ eV/q $z = 0 \pm 0.5$ mm, $E_z = 0 \pm 0.023$ eV/q correlated (i.e. ions at -0.5mm have -0.023eV/q)
 - $z = 0 \pm 0.5$ mm, $E_z = 0 \pm 0.023$ eV/q random (i.e. no correlation between energy and position).
- 500 ions used in all simulations.

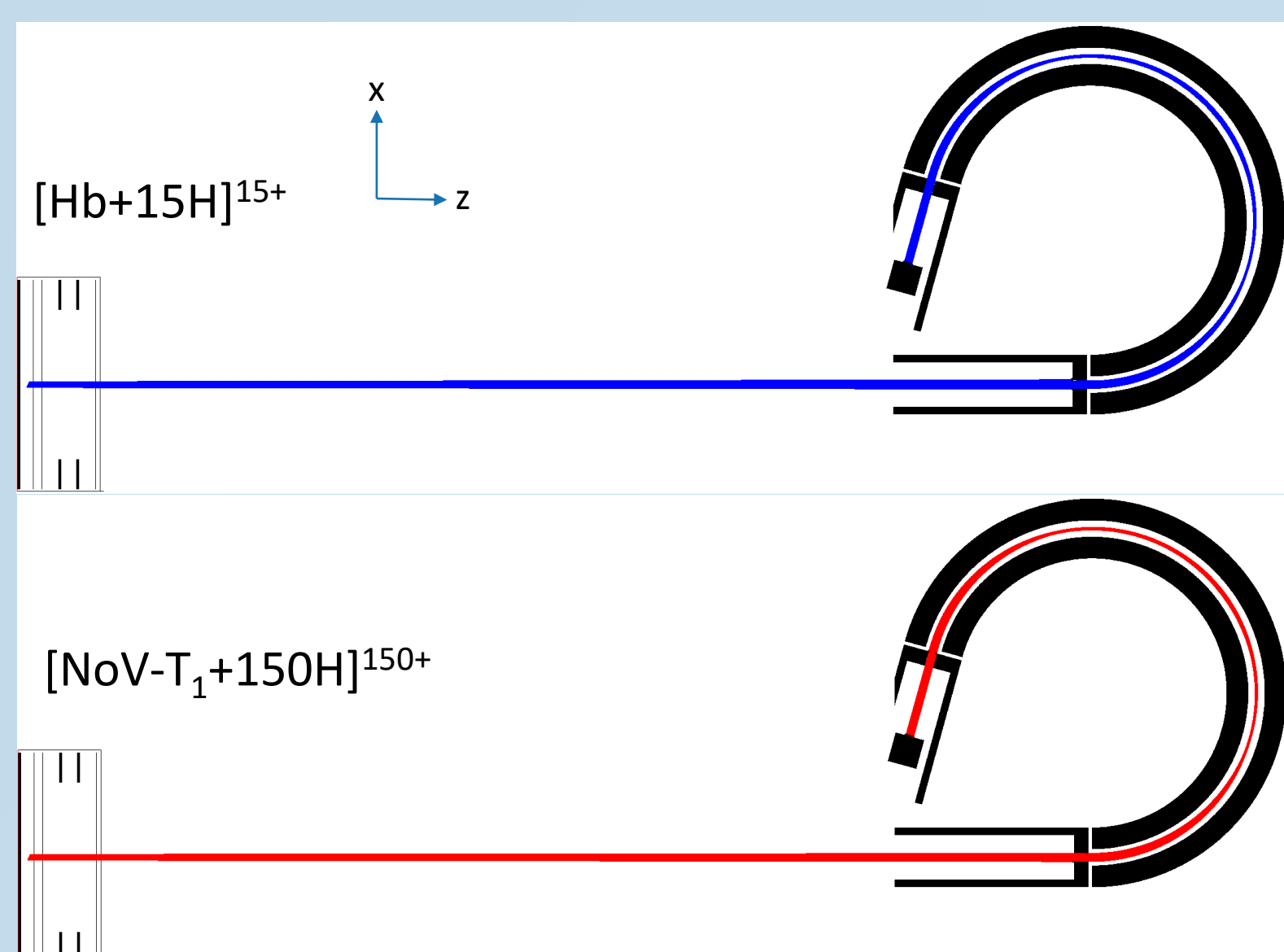

Typical ion trajectories for haemoglobin 15+ (top) and norovirus T1 capsids 150+ (bottom). Left shows the pusher stack region. Blank workbench simulated field free region. ESA shown on the right, with field free regions.

Results: Metastable Rejection

- Inherent rejection of metastables will further help resolution
- ESA is energy filter → only metastable fragments of same charge will pass.
- Possibility to scan ESA voltage to probe metastable peaks.

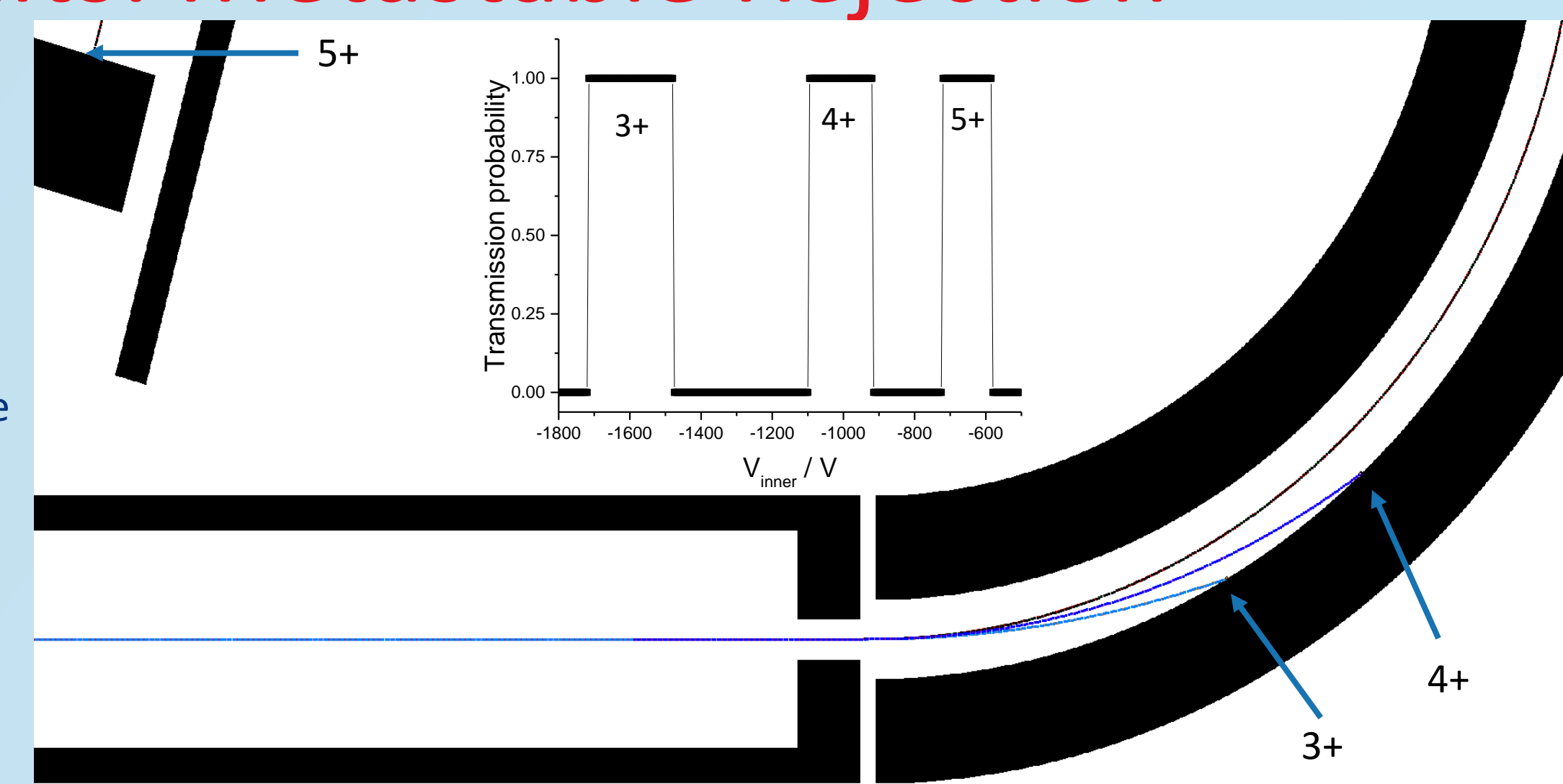

Ion trajectories for different metastable ions. Any change in E/q will prevent ions being detected. The ESA voltage can be scanned to transmit the metastable ions (inset).

Summary and Future

- Compact ESA-TOF design with a footprint less than 450 x 190 x 180 mm (lwxhx).
- Flexible design allows to alter footprint without compromising performance.
- Energy filtering of ESA removed metastable ions or allows to study them.
- Simulation of mass spectrometry performance shows high ion transmission.
- Simulated resolution > 8,000 gives excellent performance for small size.
- Prototype currently in final stages of design and ready for manufacturing.



This project has received funding from the European Union's Horizon 2020 research and innovation programme under grant agreement No 801406